


# **Class II Biological Safety Cabinets**

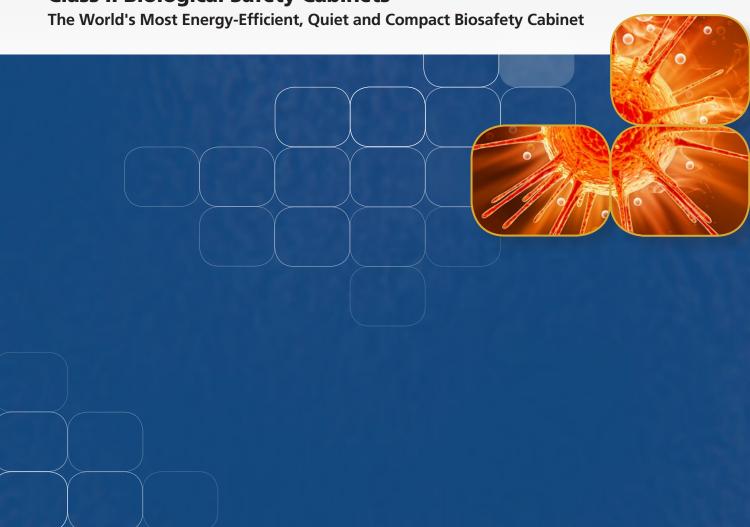



#### **RS 232 Serial Interface Port**

- Send operational information to Building Management System (BMS)
- Optional zero volt exhaust and alarm contact



B. .... AM S. HOCOL . .....

шишин

ESCO

Sentinel•GOLD







## **Sentinel™ Gold Microprocessor Controller**

- Displays all safety information on one screen
- Centered and angled down for easy reach & viewing
- Selectable Quickstart mode for fast operation



## **Angled Corner & Glass Side**

- Easy to clean
- Easy to reach service fixture and outlets
- Stainless steel side wall is available (AC2-S and AC2-D variant)



### **Divided Work Tray**

- Easy to lift and clean
- Single-piece recessed tray is available (AC2-S and AC2-D variant)



### **Raised Arm Rest**

- Helps prevent grille blocking
- Comfortable working posture







Esco Airstream® Class II has been certified by PHE / Public Health England (formerly HPA) for compliance to EN 12469

Class II Biological Safety Cabinets

### Airflow Sensor

- Monitors real-time airflow for safety
- Alerts the user if airflow is insufficient


### **Energy-Efficient DC ECM Motor**

- The most energy-efficient Class II biosafety cabinet in the world, provides 70% energy savings compared to AC motor
- Stable airflow, despite building voltage fluctuations & filter loading
- Night Setback mode to further reduce power consumption by 60%





	Biosafety Cabinet	Air Quality	Filtration	Electrical Safety
Standards Compliance	EN 12469, Europe SANS 12469, South Africa	ISO 14644.1 Class 3, Worldwide JIS B9920 Class 3, Japan JIS BS 5295, Class 3, UK	EN-1822 (H14), Europe IEST-RP-CC001.3, USA IEST-RP-CC007, USA IEST-RP-CC034.1, USA	IEC 61010-1, Worldwide EN 61010-1, Europe UL 61010-1, USA CAN / CSA-22.2, No.61010-1

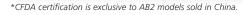





Dynamic air barrier, where inflow and forward-directed downflow air converge

ULPA-filtered air

4

- Unfiltered / potentially contaminated air
- Room air / Inflow air

### **Cabinet Filtration System**

- Ambient air is pulled through front grille to create inflow, without going through the work surface. Inflow is joined by half of the downflow, to create front air curtain that is fine-tuned to create a large performance envelope. The combined air stream travels through the back air column towards the blower.
- Approximately 1/3 of the air in the common plenum is exhausted through the ULPA filter to the room. The remaining ¾ of the air is passed through the downflow ULPA filter and into the work area as a vertical laminar flow air to create ISO Class 3 work surface and prevents cross contamination.
- Near the work surface, the downflow splits. About Half goes to the front grille, and half goes to the rear grille. A small portion enters the the side capture zones to prevent dead air corners (small blue arrows).
- The design was optimized to give large performance envelope, that provides operator and product protection at wide Inflow and Downflow variation from the Nominal point.



0.10 0.20 0.30 0.40 0.50 0.60 Downflow (m/s)

Nominal Airflow

0

- Personnel / Product Protection
- Area of Personnel / **Product Protection**
- A No Personnel / **Product Protection**
- Area of no Personnel / **Product Protection**

LCD simultaneously displays time, airflow & sash status, inflow and downflow velocities, and status remarks.

Multi Language: English, French, German, Spanish, Italian.

Diagnostics button, to easily check the cabinet operating parameters and assist servicing.

Large touchpad control buttons provide good tactile feedback. Color coded LED: green for fan; blue for FL lights and outlets; and orange for UV lamp.

Programmable UV light timer extends UV lamp life.



**Direct Mounted / GFCI** 

Outlet

Esco AC2 Airflow: OK I: 0.45 m/s

Sash: 0K D: 0.30 m/s





EO-\_









MICROPROCESSOR CONTROL SYSTEM

## **Accesories for AC2-G3 Biological Safety Cabinets**

	Glass Side Wall		AC2-2E8 <b>2010718</b>	AC2-3E8 <b>2010658</b>	AC2-4E8 <b>2010621</b>	AC2-5E8 <b>2010656</b>	AC2-6E8 <b>2010657</b>	
		230 VAC,			AC2-4G8 <b>2010734</b>		AC2-6G8 <b>2010743</b>	
		50/60 Hz	AC2-2S8 <b>2010767</b>	AC2-3S8 <b>2010721</b>	AC2-4S8 <b>2010711</b>	AC2-5S8 <b>2010725</b>	AC2-6S8 <b>2010722</b>	
Cabinet	Stainless Steel Side Wall				AC2-4D8 <b>2010733</b>		AC2-6D8 <b>2010742</b>	
	Glass Side Wall	115 VAC, 50/60 Hz	AC2-2E9 <b>2010777</b>	AC2-3E9 <b>2010779</b>	AC2-4E9 <b>2010697</b>	AC2-5E9 <b>2010784</b>	AC2-6E9 <b>2010787</b>	
	Stainless Steel Side Wall		AC2-2S9 <b>2010790</b>	AC2-3S9 <b>2010792</b>	AC2-4S9 <b>2010744</b>	AC2-5S9 <b>2010797</b>	AC2-6S9 <b>2010800</b>	
	Anti-blowback Valve 10"		ABBV-10P 5170352					
Exhaust Ducting	Tri-safe Exhaust Collar with Alarm		N/A		TEM-4 N/A 2010606		/A	
	Thimble Exhaust Collar		ECO-AC22 5170520	ECO-AC23 5170521	ECO-AC24 5170522	ECO-AC25 5170523	ECO-AC26 5170524	
	UV Lamp		UV-15A-L 5170251		UV-30A-L 5170255			
	IV Bar		IV-605 5170498	IV-910 5170499	IV-1215 5170231	IV-1520 5170500	IV-1825 5170501	
Work Zone	Multiple Piece Tray Option (AC2-5 / AC2-D)		SDT-AC2-2E 5020643	SDT-AC2-3E 5020635	SDT-AC2-4E 5020606	SDT-AC2-5E 5020640	SDT-AC2-6E 5020592	
	Single Piece Tray Option (AC2-E / AC2-G)		SGT-AC2-2S 5020696	SGT-AC2-3S 5020648	SGT-AC2-4S 5020627	SGT-AC2-5S 5020651	SGT-AC2-6S 5020645	
	Pre-filter		PF-40 5090060	PF-41 5090061	PF-42 5090062	PF-43 5090063	PF-44 5090064	
Electrical	ical Direct Mounted / GFCI				FO-			



ABBV-10P



TEM-4



ECO-AC2-G3



Service Fixtures	EU SF-Gas-20 mm		S	F-1G20 517041	0				
	EU SF-Vacuum-20 mm	SF-1V20 5170457							
	EU SF-Air-20 mm	SF-1A20 5170502							
	EU SF-Nitrogen-20 mm		SF-1N20 5170503						
	EU SF-Water-20 mm	SF-1W20 5170458							
	EU SF-Universal-22 mm	SF-2U22 5170504							
	Fixed Stand with Leveling Feet, 28" height	SAL-2A0 Gen 2 5130169	SAL-3A0 Gen 2 5130170	SAL-4A0 Gen 2 5130134	SAL-5A0 Gen 2 5130171	SAL-6A0 Gen 2 5130172			
	Fixed Stand with Leveling Feet, 34" height	SAL-2B0 Gen 2 5130173	SAL-3B0 Gen 2 5130174	SAL-4B0 Gen 2 5130175	SAL-5B0 Gen 2 5130176	SAL-6B0 Gen 2 5130177			
	Fixed Stand with Caster Wheels, 28" height	SPC-2A0 Gen 2 5130161	SPC-3A0 Gen 2 5130155	SPC-4A0 Gen 2 5130152	SPC-5A0 Gen 2 5130162	SPC-6A0 Gen 2 5130154			
Support Stands, Ships Flat	Fixed Stand with Caster Wheels, 34" height	SPC-2B0 Gen 2 5130164	SPC-3B0 Gen 2 5130165	SPC-4B0 Gen 2 5130166	SPC-5B0 Gen 2 5130167	SPC-6B0 Gen 2 5130168			
Snips riat	Telescopic Stand with Leveling Feet, 1" adjustment	STL-2A0 5130092	STL-3A0 5130050	STL-4A0 5130051	STL-5A0 5130052	STL-6A0 5130053			
	Telescopic Stand with Caster Wheels, 1" adjustment	STC-2A0 5130135	STC-3A0 5130055	STC-4A0 5130056	STC-5A0 5130057	STC-6A0 5130058			
	Motorized Height Stand with Caster Wheels, 39.5" height		SPM-3A2 5130093	SPM-4A2 5130047	SPM-5A2 5130100	SPM-6A2 5131141			
Misc	IQ/OQ Protocol	9010179							



PF-\_



ABBV-10P



SF-





SPC-\_A0 Gen2

15

5

### **AC2 Biological Safety Cabinet Engineering Drawing**

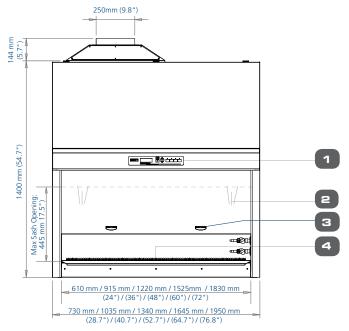


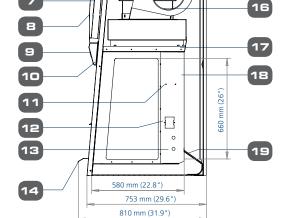

- 1. Esco Sentinel™ Gold Microprocessor Control System
- 2. Tempered Glass Sliding Sash Window
- 3. Ergonomic Sash Handle
- 4. Stainless Steel Work Tray (available in single and multi pieces)
- 5. Exhaust Collar (optional)
- 6. Exhaust ULPA / H14 Filter

- 7. Electrical Panel
- 8. Curved Front Panel
- 9. Angled down LCD and Control Panel
- 10. T5 Fluorescent Lamps
- 11. IV Bar Retrofit Kit Provision
- 12. Electrical outlet Retrofit Kit Provision (1 on each side)
- 13. Service Fixture Retrofit Kit Provision (2 on each side)
- 14. Ergonomic Dual Posture Stainless Steel Arm Rest
- 15. DC ECM Blower
- 16. Second Blower, for exhaust (AC2-D & AC2-G)
- 17. Downflow ULPA / H14 Filter
- 18. UV Lamp (optional)
- 19. Removable Paper Catch (with optional pre-filter)

## **Optional Exhaust Collar Positions for Thimble-Ducting for AC2 Models**

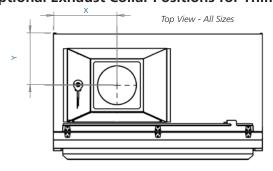

Size	2	3	4	5	6	ft
	0.6	0.9	1.2	1.5	1.8	m
х	233	331	408	560	560	
Y	334	334	334	334	326	mm
х	9.2	13	16	22	22	inches
Y	13.1	13.1	13.1	13.1	12.8	inches

		TECHNIC	AL SPECIFICATIO	NS					
Glass Side: 230 VAC, 50/60 Hz		AC2-2E8 2010718	AC2-3E8 2010658	AC2-4E8 2010621	AC2-5E8 2010656	AC2-6E8 2010657			
Glass Side: 115 VAC, 50/60 Hz		AC2-2E9 2010777	AC2-3E9 2010779	AC2-4E9 2010697	AC2-5E9 2010784	AC2-6E9 2010787			
Stainless Steel Side: 230 VAC, 50/60 Hz		AC2-258 2010767	AC2-358 2010721	AC2-458 2010711	AC2-5S8 2010725	AC2-658 2010722			
Stainless Steel Side: 11	15 VAC, 50/60 Hz	AC2-2S9 2010790	AC2-359 2010792	AC2-4S9 2010744	AC2-5S9 2010797	AC2-659 2010800			
Nominal Size		2 ft (0.6 meter)	3 ft (0.9 meter)	4 ft (1.2 meter)	5 ft (1.5 meter)	6 ft (1.8 meter)			
	Width	730 mm (28.8")	1035 mm (40.8")	1340 mm (52.8")	1645 mm (64.8")	1950 mm (76.8")			
External Dimensions	Depth without Arm Rest		<u> </u>	753 mm (29.5")		<u>I</u>			
(W x D x H)	Depth with Arm Rest	810 mm (32.0")							
	Height	1400 mm (54.8")							
	Width	610 mm (24.0")	915 mm (36.0")	1220 mm (48.0")	1525 mm (60.0")	1830 mm (72.0")			
Gross Internal Dimensions	Depth			580 mm (22.8")		<u> </u>			
(W x D x H)	Height			660 mm (26.0")					
Usable Work Area		0.27 m <sup>2</sup> (2.9 sq.ft.)	0.42 m <sup>2</sup> (4.5 sq.ft.)	0.56 m <sup>2</sup> (6.1 sq.ft.)	0.71 m² (7.63 sq.ft.)	0.86 m² (9.2 sq.ft.)			
Tested Opening				175 mm (7")					
Work Opening				190 mm (7.5")					
Maximum Opening		475 mm (18.7")							
Inflow		0.45 m/s (90 fpm)							
Average Airflow Velocity	Downflow	0.30 m/s (60 fpm)							
	Inflow	173 cmh (102 cfm)	259 cmh (152 cfm)	346 cmh (204 cfm)	432 cmh (254 cfm)	519cmh (305 cfm)			
	Downflow	369 cmh (217 cfm)	553 cmh (325 cfm)	738 cmh (434 cfm)	922 cmh (543 cfm)	1107 cmh (657 cfm)			
	Exhaust	173 cmh (102 cfm)	259 cmh (152 cfm)	346 cmh (204 cfm)	432 cmh (254 cfm)	519cmh (305 cfm)			
Airflow Volume	Required Exhaust with Optional Thimble Exhaust Collar	260 m³/h (153 cfm)	320 m³/h (189 cfm)	538 m³/h (317 cfm)	615 m³/h (362 cfm)	823 m³/h (485 cfm)			
	Static Pressure for Optional Thimble Exhaust Collar	28 Pa / 0.11 in H <sub>2</sub> O	29 Pa / 0.11 in H <sub>2</sub> O	31 Pa / 0.12 in H <sub>2</sub> O	35 Pa / 0.14 in H <sub>2</sub> O	47 Pa / 0.18 in H <sub>2</sub> O			
III DA Eiltor Tunical Effic	ioner	>99.999% at 0.1 to 0.3 micron, ULPA as per IEST-RP-CC001.3 USA							
ULPA Filter Typical Effic	iency	>99.999% at MPPS, H14 as per EN 1822 EU							
Sound Emission*	NSF / ANSI 49	56.3	56.6	58.7	58.2	59.4			
Souria Emission"	EN 12469	51.0	52.0	53.5	53.6	55.7			
Fluorescent Lamp Inten	sity (Lux)	859	1279	1404	1227	1384			
Fluorescent Lamp Inten	sity (foot-candles)	80	119	130	114	129			
	Main body	1.2 mm (0.05") 18 gauge electro-galvanized steel with white oven-baked epoxy-polyester Isocide™ antimicrobial powder-coated finish							
Cabinet Construction	Work Zone		1.5 mm (0.06") 16 g	gauge stainless steel, type	e 304, with 4B finish				
	Side Walls (E Series)		UV-absorbing tempere	ed glass, 5 mm (0.2"), col	orless and transparent				
	Side Walls (S Series)	1.5 mm (0.06") 16 gauge stainless steel, type 304, with 4B finish							
Electrical	Cabinet Full Load Amps (FLA)	1.8	3.5	3.7	4.3	5.5			
Electrical	Heat Load (BTU/Hr)	324	447	580	717	966			
Nominal Power Consum	nption (W)	95	131	160	210	283			
Net Weight**		116 Kg (256 lbs)	173 Kg (381 lbs)	230 Kg (507 lbs)	288 Kg (635 lbs)	346 Kg (763 lbs)			
Shipping Weight**		143 Kg (315 lbs)	214 Kg (472 lbs)	285 Kg (628 lbs)	356 Kg (785 lbs)	428 Kg (944 lbs)			
Shipping Dimensions Maximum (W x D x H)*	*	850 x 820 x 1760 mm (33.5" x 32.3" x 69.3")	1120 x 820 x 1760 mm (44.1" x 32.3" x 69.3")	1450 x 820 x 1760 mm (57.1" x 32.3" x 69.3")	1720 x 820 x 1760 mm (67.7" x 32.3" x 69.3")	2050 x 820 x 1760 mm (80.7" x 32.3" x 69.3")			
Shipping Volume, Maxi	imum**	1.23 m³ (43.2 ft³)	1.62 m³ (57.2 ft³)	2.09 m <sup>3</sup> (73.8 ft <sup>3</sup> )	2.48 m <sup>3</sup> (87.6 ft <sup>3</sup> )	2.96 m³ (104.5 ft³)			

<sup>\*</sup>Noise reading in open field condition / **anechoic** chamber. Noise reading in **normal room varies** by room size, layout, and background noise, but may reach roughly 3-4 dBA above these values. \*\*Cabinet only, excludes optional stand.

TECHNICAL SPECIFICATIONS					
Glass Side: 230 VAC, 50/60 Hz		AC2-4G8 2010734	AC2-6G8 2010743		
Stainless Steel Side: 23	0 VAC, 50/60 Hz	AC2-4D8 2010733	AC2-6D8 2010742		
Nominal Size		4 ft (1.2 meter)	6 ft (1.8 meter)		
	Width	1340 mm (52.8")	1950 mm (76.8")		
External Dimensions	Depth without Arm Rest	753 mm	1 (29.5")		
(W x D x H)	Depth with Arm Rest	810 mm	1 (32.0")		
	Height	1400 mm (54.8")			
Gross Internal	Width	1220 mm (48")	1830 mm (72")		
Dimensions	Depth	580 mm	1 (22.8")		
(W x D x H)	Height	660 mm (26")			
Usable Work Area		0.56 m² (6.1 sq.ft.)	0.86 m² (9.0 sq.ft.)		
Tested Opening		175 m	m (7")		
Working Opening		190 mr	n (7.5")		
Average Airflow	Inflow	0.45 m/s	(90 fpm)		
Velocity	Downflow	0.30 m/s	(60 fpm)		
	Inflow	346 cmh (588 cfm)	519 cmh (881 cfm)		
	Downflow	738 cmh (1254 cfm)	1107 cmh (1880 cfm)		
	Exhaust	346 cmh (588 cfm)	519 cmh (881 cfm)		
Airflow Volume	Required Exhaust with Optional Thimble Exhaust Collar	538 m³/ h (317 cfm)	823 m³/ h (485 cfm)		
	Static Pressure for Optional Thimble Exhaust Collar	31 Pa / 0.12 in H <sub>2</sub> O	47 Pa / 0.18 in H <sub>2</sub> O		
ULPA Filter Typical Effici	encv	>99.999% at 0.1 to 0.3 micron, l	JLPA as per IEST-RP-CC001.3 USA		
OLI ATTICCI Typical Effici	-	>99.999% at MPPS, H14 as per EN 1822 EU			
Sound Emission*	NSF / ANSI 49	61.3 dBA	62.5 dBA		
Journa Emission	EN 12469	58.3 dBA	59.5 dBA		
Fluorescent Lamp Inten	sity (Lux)	1400			
Fluorescent Lamp Inten	sity (foot-candles)	13	30		
	Main body	1.2 mm (0.05") 18 gauge electro-galvanized steel with white oven-baked epoxy-polyester Isocide™ antimicrobial powder-coated f			
Cabinet Construction	Work Zone	1.5 mm (0.06") 16 gauge stainle	ess steel, type 304, with 4B finish		
	Side Walls (G-Series)	UV-absorbing tempered glass, 5 m	m (0.2"), colorless and transparent		
	Side Walls (D-Series)	1.5 mm (0.06") 16 gauge stainle	ess steel, type 304, with 4B finish		
Florendes	Cabinet Full Load Amps (FLA)	9.4 A	12.6 A		
Electrical	Heat Load (BTU / Hr)	905	1230		
Nominal Power Consum	nption	265 W	360 W		
Net Weight**		240 Kg (529 lbs)	366 Kg (807 lbs)		
Shipping Weight**		295 Kg (650 lbs)	448 Kg (988 lbs)		
Shipping Dimensions, N	flaximum (W x D x H)**	1450 x 820 x 1760 mm (57.1" x 32.3" x 69.3")	2050 x 820 x 1760 mm (80.7" x 32.3" x 69.3")		
Shipping Volume, Maxi	mum**	2.09 m³ (73.8 ft³)	2.96 m³ (104.5 ft³)		

\*Noise reading in open field condition / anechoic chamber. Noise reading in normal room varies by room size, layout, and background noise, but may reach roughly 3-4 dBA above these values. \*\*Cabinet only, excludes optional stand.

First Airstream® Offers the Most Complete Class II Cabinet Range


1.5 m (5'), 1.8 m (6')

Available

Airstream Product	E-Series G-Series		S-Series	D-Series	
		One-piece stainless steel with coved corners for cleanability. Side captures and negative pressure side walls optimize containment.			
Work Tray	Multi-piece, Autoclavable		Single-piece stainless steel, spill retaining		
Fan System	Single blower for inflow and downflow. Energy-efficient and cost-effective	Dual blowers for inflow and downflow. Redundant system provides protection in case of fan failure.	Single blower for inflow and downflow. Energy-efficient and cost-effective.	Dual blowers for inflow and downflow. Redundant system provides protection in case of fan failure.	
Exhaust Filter	Filter with > 99.999% efficiency > 100x better protection than single filter system  Size 0.6 m (2'), 0.9 m (3'), 1.2 m (4'), 1.2 m (4') 1.8m (6')		Single ULPA/H14 Filter >99.999% Efficient, Cost Effective	Dual ULPA/H14 Filters, >100.000x Better Protection than Single Filter System	
Size			0.6 m (2'),0.9 m (3'), 1.2 m (4'),	1.2 m (4'), 1.8 m (6')	


1.5 m (5'), 1.8 m (6')

### **AC2-D Airflow Diagram**

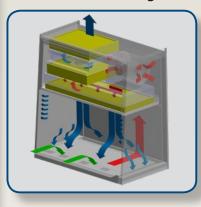

- ULPA-filtered air
- Unfiltered / potentially contaminated air
- Room air / Inflow air

### **AC2-G Airflow Diagram**

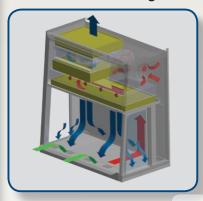

- ULPA-filtered air
- Unfiltered / potentially contaminated air
- Room air / Inflow air

## Comprehensive **Performance Testing At Esco**

Every Airstream® AC2 model manufactured by Esco is individually tested, documented by serial number and validated with the following test methods:

- Inflow and downflow velocity
- PAO aerosol challenge for filter integrity
- Airflow pattern visualization
- Electrical safety to IEC61010-1
- Additional KI-Discus containment and microbiological testing are performed on statistical sampling basis.





#### **ESCO GLOBAL NETWORK**

42 LOCATIONS IN 21 COUNTRIES ALL OVER THE WORLD



Biological Safety Cabinets

CO<sub>2</sub> Incubators

Compounding Pharmacy Equipment
Containment / Pharma Products

Ductless Fume Hoods

Lab Animal Research Products

Laboratory Centrifuges

Laboratory Fume Hoods

Laboratory Ovens and Incubators

Laboratory Shakers

Laminar Flow Clean Benches

PCR Cabinets

PCR Thermal Cyclers

Powder Weighing Balance Enclosures

Ultra-low Temperature Freezers

The Esco Group of Companies is a global life sciences tools provider with sales in over 100 countries. The group is active in lab equipment, pharma equipment and medical devices. Manufacturing facilities are located in Asia and Europe. R&D is conducted worldwide spanning the US, Europe and Asia. Sales, service and marketing subsidiaries are located in 12 major markets including the US, UK, Singapore, Japan, China and India. Regional distribution centers are located in the US, UK, and Singapore.

Life Science • Chemical Research • Assisted Reproductive Technology (ART) • Pharmaceutical Equipment • General Equipment



### WORLD CLASS, WORLDWIDE.

Esco Micro Pte. Ltd. • 21 Changi South Street 1 • Singapore 486 777 Tel +65 6542 0833 • Fax +65 6542 6920 • mail@escoglobal.com www.escoglobal.com

Esco Technologies, Inc. • 903 Sheehy Drive, Suite F, Horsham, PA 19044, USA Tel 215-441-9661 • Fax 484-698-7757 eti.admin@escoglobal.com • www.escolifesciences.us

Esco Global Offices: Bangladesh | China | Denmark | Germany | Hong Kong | Indonesia | Italy | Lithuania | Malaysia | Myanmar | Philippines | Russia | Singapore | South Africa | South Korea | Taiwan | Thailand | UAE | UK | USA | Vietnam





